# Package: musicNMR (via r-universe)

September 7, 2024

Version 1.0

Date 2024-01-09

Maintainer Stefano Cacciatore <tkcaccia@gmail.com>

Title Conversion of Nuclear Magnetic Resonance Spectra in Audio Files

**Description** A collection of functions for converting and visualization the free induction decay of mono dimensional nuclear magnetic resonance (NMR) spectra into an audio file. It facilitates the conversion of Bruker datasets in files WAV. The sound of NMR signals could provide an alternative to the current representation of the individual metabolic fingerprint and supply equally significant information. The package includes also NMR spectra of the urine samples provided by four healthy donors. Based on Cacciatore S, Saccenti E, Piccioli M. Hypothesis: the sound of the individual metabolic phenotype? Acoustic detection of NMR experiments. OMICS. 2015;19(3):147-56. <doi:10.1089/omi.2014.0131>.

**Depends** R ( $\geq$  2.10.0), seewave

SuggestsNote No suggestions

License GPL (>= 2)

NeedsCompilation no

Date/Publication 2024-01-10 13:53:19 UTC

Author Stefano Cacciatore [aut, trl, cre] (<https://orcid.org/0000-0001-7052-7156>)

Repository https://tkcaccia.r-universe.dev

RemoteUrl https://github.com/cran/musicNMR

RemoteRef HEAD

RemoteSha faa577db7a2d3c9f45fd51d7627b97b7716bfcef

# Contents

AU..... 2

Α	Ψ	 •													•						3	1
m	usicMatrix	 •													•						4	
m	usicNMR	 •													•						5	
pl	otFID	 •													•						6	)
																					7	

# Index

AU

Collection of Free Induction Decay of Urine Spectra of the Donor AU

#### Description

This dataset consists of eight Free Induction Decay (FID) from eight urine samples. All urine were collected from the donor "AU". The urine are divided in two group; in each group urine samples were collected in a time period of a week. The collection of the two groups is separated by a time course of two years.

# Usage

data(AU)

#### Value

A list with the following elements:

sample1A	a matrix object. FID of the sample 1 of group A.
sample2A	a matrix object. FID of the sample 2 of group A.
sample3A	a matrix object. FID of the sample 3 of group A.
sample4A	a matrix object. FID of the sample 4 of group A.
sample1B	a matrix object. FID of the sample 1 of group B.
sample2B	a matrix object. FID of the sample 2 of group B.
sample3B	a matrix object. FID of the sample 3 of group B.
sample4b	a matrix object. FID of the sample 4 of group B.

# References

Assfalg M, Bertini I, Colangiuli D, *et al.* Evidence of different metabolic phenotypes in humans. *Proc Natl Acad Sci U S A* 2008;105(5):1420-4.

### Examples

```
data(AU)
plotFID(AU$sample1B)
plotFID(AU$sample1A,ADD=TRUE,col=2)
```

# Description

This dataset consists of eight Free Induction Decay (FID) from eight urine samples. All urine were collected from the donor "AW". The urine are divided in two group; in each group urine samples were collected in a time period of a week. The collection of the two groups is separated by a time course of two years.

#### Usage

data(AW)

# Value

A list with the following elements:

sample1A	a matrix object. FID of the sample 1 of group A.
sample2A	a matrix object. FID of the sample 2 of group A.
sample3A	a matrix object. FID of the sample 3 of group A.
sample4A	a matrix object. FID of the sample 4 of group A.
sample1B	a matrix object. FID of the sample 1 of group B.
sample2B	a matrix object. FID of the sample 2 of group B.
sample3B	a matrix object. FID of the sample 3 of group B.
sample4b	a matrix object. FID of the sample 4 of group B.

# References

Assfalg M, Bertini I, Colangiuli D, *et al.* Evidence of different metabolic phenotypes in humans. *Proc Natl Acad Sci U S A* 2008;105(5):1420-4.

# Examples

```
data(AW)
plotFID(AW$sample1B)
plotFID(AW$sample1A,ADD=TRUE,col=2)
```

# AW

musicMatrix

# Description

This function save a matrix in audio file as \*.wav

# Usage

musicMatrix(ma,destination)

#### Arguments

ma	a matrix. The time in second unit is in the firt column. The intensity of the FID
	is in the second column.
destination	Name of the .wav file

#### Details

This function uses - three functions from the package **tuneR**: Wave, normalize and writeWave; - one function from the package **seewave**: savewav.

# Value

No return value.

# Author(s)

Stefano Cacciatore

#### References

Cacciatore S, Saccenti E, Piccioli M. Hypothesis: the sound of the individual metabolic phenotype? Acoustic detection of NMR experiments. OMICS. 2015 Mar;19(3):147-56.

# Examples

```
data(AU)
musicMatrix(AU$sample1A,"audio.wav")
```

# The function unlink deletes the file unlink("audio.wav") musicNMR

#### Description

This function converts monodimensional Nuclear Magnetic Resonance spectrum in audio file. The spectrum is imported as Bruker format and the audio file is saved as \*.wav

# Usage

musicNMR(source,destination)

#### Arguments

source	The folder's address where the Free Induction Decay is located
destination	Name of the .wav file

#### Details

This function uses - three functions from the package **tuneR**: Wave, normalize and writeWave; - one function from the package **seewave**: savewav.

#### Value

No return value.

#### Author(s)

Stefano Cacciatore

#### References

Cacciatore S, Saccenti E, Piccioli M. Hypothesis: the sound of the individual metabolic phenotype? Acoustic detection of NMR experiments. OMICS. 2015 Mar;19(3):147-56.

plotFID

# Description

A function for ploting Free Induction Decay (FID).

# Usage

plotFID(x,ADD=FALSE, ...)

# Arguments

x	a matrix. The time in second unit is in the firt column. The intensity of the FID is in the second column.
ADD	For a new plot ADD = FALSE. To overimpose a different FID on an existing plot ADD = TRUE.
	Arguments to be passed to the plot function.

# Value

No return value.

# Author(s)

Stefano Cacciatore

#### References

Cacciatore S, Saccenti E, Piccioli M. Hypothesis: the sound of the individual metabolic phenotype? Acoustic detection of NMR experiments. OMICS. 2015 Mar;19(3):147-56.

# Examples

```
data(AU)
plotFID(AU$sample1B)
plotFID(AU$sample1A,ADD=TRUE,col=2)
```

# Index

\* datasets AU, 2 AW, 3 AU, 2 AW, 3 musicMatrix, 4 musicNMR, 5 normalize, 4, 5 plot, 6 plotFID, 6 savewav, 4, 5 Wave, 4, 5 writeWave, 4, 5